MICROTECH SX93 Stereo Microscope Operation Manual

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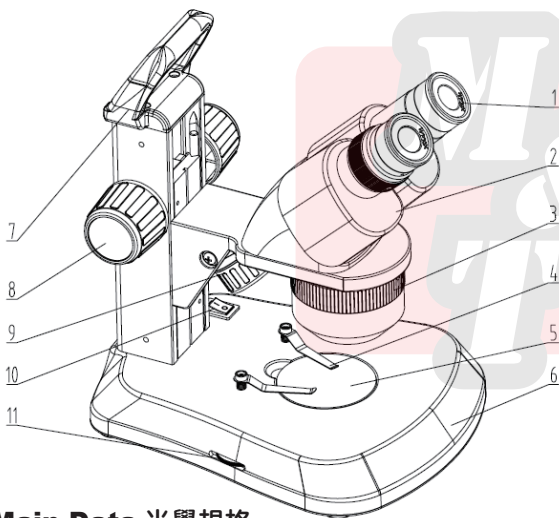
Please read carefully before using the microscope

1. Applications 應用

Widely used in electronical industry, assembly and inspection of precision instrument, dissection and study of propagation, also applied in field of school, carving, geology, archeology etc
廣泛的應用於電子工業、裝配維修精密儀器、研究檢驗、生物解剖、學術、地質、考古、美術雕刻、珠寶鑲埋、寶石鑑定、顯微鑑識等，用途極為廣泛。

2. Name of the parts and accessories 部位說明

1. Eyepiece 目鏡
2. Prism Box 稜鏡室
3. Objective cover 物鏡
4. Clips 固定夾
5. Working stage 工作台附底光
6. Base底座
7. Handle 手提把
8. Focusing Knob 調焦鈕
9. Incident Illumination 投射照明
10. Power Switch 電源開關
11. Dimmer 調光鈕

3. Main Data 光學規格

1) Optical Data (mm)

Eyepiece 目鏡/倍數		WF 5X	WF 10X	WF 15X	WF 20X	
Objective 物鏡倍率	1X 標配	Field of View視野mm	22	20	15	10
		Working Distance工作距離mm	57	57	57	57
	2X 標配	Field of View視野mm	12	10	7.5	6.5
		Working Distance工作距離mm	80	80	80	80
	3X	Field of View視野mm	8	6.7	5	4.3
		Working Distance工作距離mm	61	61	61	61
	4X 標配	Field of View視野mm	6	5	3.7	3.2
		Working Distance工作距離mm	57	57	57	57

(Note: Objective:1X/2X/3X or 1X/2X/4X. Optional eyepiece WF5X,WF15X,WF20X)

5. LED lamp and battery replacement 燈泡與電池的更換

1) Replace the incident bulb 更換投射燈泡

Firstly remove objective cover via taking off 4 screws. Secondly take off the lampshade, replace the burned bulb with new bulb by input the feet to the porcelain socket fully. Then set the lampshade on again.

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2) Replace the transmitted bulb 更換穿透光燈泡

Firstly loosen the screw, then take off frosted glass and filter, take out the burned bulb and replace with new bulb by input the feet on the porcelain socket fully, at last set back the frosted glass and screw.

Note: Usually longer pins means anode.

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3) Replace battery 更換電池

Open battery box under base of microscope, take out old batteries and replace new batteries. Please put out

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6. Maintaining and storage 維修及存放

- 1) Microscope is a precision instrument, you should place carefully, avoid impact during transportation.
- 2) Put it in a dry and clean place, avoid high temperature and shock.
- 3) Do not touch the lens directly by your finger.
- 4) Keep the optical surface clean, dust on the surface you can wipe off by hair drier.
- 5) Do not use organic things to wipe the surface of microscope, especially plastic surface, please clean with neutral scour.
- 6) All the optical and mechanism part are adjusted and could not take apart by yourself.
- 7) Add grease on moving part regularly.
- 8) Put the microscope in a cool and dry place and cover it with dust cover when not be used for a long time.

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3) Structure data 結構規格

Eyepiece目鏡：WF10X(標配) Objective物鏡倍率(三組)：1X/2X/4X

45° inclined binocular head, left ocular-tube is with diopter adjustment ± 5 dp, the interpupillary distance is between 54~76mm

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2) Power control 電源控制

Turn on the switch, adjust the brightness of Incident and Transmitted light by the two dimmers.

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3) Selecting of working stage 工作台的選擇

a) Normally frosted glass stage equipped with microscope when packing, locked by screw. You can use transmit light to observe transparent object. 使用毛玻璃時可以利用穿透光觀察透明物體。

b) Black and white working stage is packed with the microscope accessories for choice. When use, please loosen the screw and take off the glass stage. Normally use the white side. For white or other transparent object, you need to use black side to improve the contrast with Incident light. 使用黑白不透光工作台，通常使用白色面，若觀察白色或透明物體則使用黑色面，可以有效改善對比度與光的反射。

4) Usage of eye's cover 眼罩的使用

The eye's covers are packed with the microscope as accessory. Please put them on the eyepiece when you need. 橡皮製眼罩是顯微鏡附件，適合未戴眼鏡觀察者，建議安裝在目鏡上使用，可有效保持目鏡與眼球的距離。

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Place the specimen on the centre of the working stage, rotate magnification knob (or objective cover) to a high power, then turn the knob slowly till you can see clearly image on right eyepiece; Then observe through left eyepiece, adjust diopter if not clear, then turn around the left and right prism house till the interpupillary distance are suitable. If need to change the magnification, just turn around magnification knob (or the objective cover). If necessary, you can adjust head's height by locking knob.

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6) Battery and adapter 電池和變壓器

This model use battery and AC adapter for power supply. 使用電池和交流電適配器電源。

When use battery, please open battery box and put 3pcs batteries in it. Then close the battery box.

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M&T OPTICS CO.,LTD